Deoxyribonuclease I (DNase I) **RNase-free**

Nonspecific deoxyribonuclease that degrades both double-stranded and single-stranded DNA endonucleolytically releasing 5'-phosphorylated di-, tri-, and oligonucleotide products.²

Cat. No.	Size
E1345-01	1 000 units
E1345-02	5 000 units

Applications:

- Preparation of DNA-free RNA (degradation of contaminating DNA after RNA isolation).³
- Preparation of DNA-free RNA prior RT-PCR and RT-qPCR.⁴
- Removal of template DNA following in vitro transcription.
- Studies of DNA-protein interactions (footprinting).
- DNA labeling by nick-translation.
- Production of random fragments (generation of libraries).⁵

of enzyme required to completely degrade 1 µg of plasmid DNA in 10 min at 37°C.

One functional Dnase I unit is unit.1

Unit Definition: One unit is the amount Enzyme activity: DNase I requires Ca²⁺ and Mg²⁺ for hydrolyzing double-stranded DNA. In the presence of Mg^{2+,} DNase I cleaves each strand of double-stranded DNA independently in a statistically random fashion (recommended Reaction Buffer I). In the presence of Mn^{2+} , the enzyme cleaves both DNA strands at approximately the same site, approximately equivalent to 0.3 Kunitz producing DNA fragments with blunt-ends or with overhang termini of only one or two nucleotide (recommended Reaction Buffer II).⁵

Storage Conditions: Store at -20°C

Quality Control: Functionally tested for digesting of template - plasmid DNA. The absence of RNase confirmed by appropriate quality test utilizing spectrophotometry of RNA assays sample concentration before and after incubation with an excess of enzyme.

Note 1: This DNase solution does not contain an RNase inhibitor. Please handle with care to avoid contaminating it with RNase.

Note 2: DNase I is sensitive to physical denaturation. Therefore, do not vortex solutions containing DNase I, mix by gently flipping the tube or pipetting.

10 x Reaction Buffer I:

100 mM Tris-HCl, 25 mM MgCl₂, 100 mM CaCl₂, pH 7.4 @ 25°C.

10 x Reaction Buffer II:

100 mM Tris-HCl, 100 mM CaCl₂, pH 7.4 @ 25°C. Supplemented with 100 mM MnCl_{2.}

Inactivation: Inactivated by heating at 65°C for 10 min in the presence of EDTA or EGTA.

Inhibitors: metal chelators (EGTA, EDTA), transition metals, SDS, reducing agents (DTT, β -mercaptoethanol).

References:

- Kunitz, M (1950) J. Gen Physiol 33: 349-362. 1.
- 2. Vanecko, S and Laskowski, M (1961). J Biol Chem 236: 3312-3316.
- Although it is not required for most applications please see the 3. additional DNase I digestion conditions in Manual for GeneMATRIX UNIVERSAL DNA/RNA/Protein Purification Kit (E3597) and GeneMATRIX UNIVERSAL RNA Purification Kit (E3598).
- 4. Sanyal, A., et al., An effective method of completely removing contaminating genomic DNA from an RNA sample to be used for PCR. Mol. Biotechnol., 8, 135-137, (1997).
- 5. Sambrook, J., Fritsch, E.F. and Maniatis, T. (1989) Molecular Cloning: A Laboratory Manual 2nd ed., Cold Spring Harbor Laboratory, Cold Spring Harbor, New York.

This product is developed, designed and sold exclusively for research purposes and in vitro use only.

EURx Ltd. 80-297 Gdansk Poland ul. Przyrodnikow 3, NIP 957-07-05-191, KRS 0000202039, www.eurx.com.pl orders: orders@eurx.com.pl tel. +48 58 524 06 97, fax +48 58 341 74 23